


LabChart 7 Feature Manager allows users to browse and download any useful ADInstruments software in one dialog. Softwares are neatly categorized according to applications to improve ease of use. Software updates can be manually or automatically performed to alert users about new versions of LabChart, LabChart Modules and LabChart Extensions.

**Note:** Only registered PowerLab users can download LabChart products and software updates.

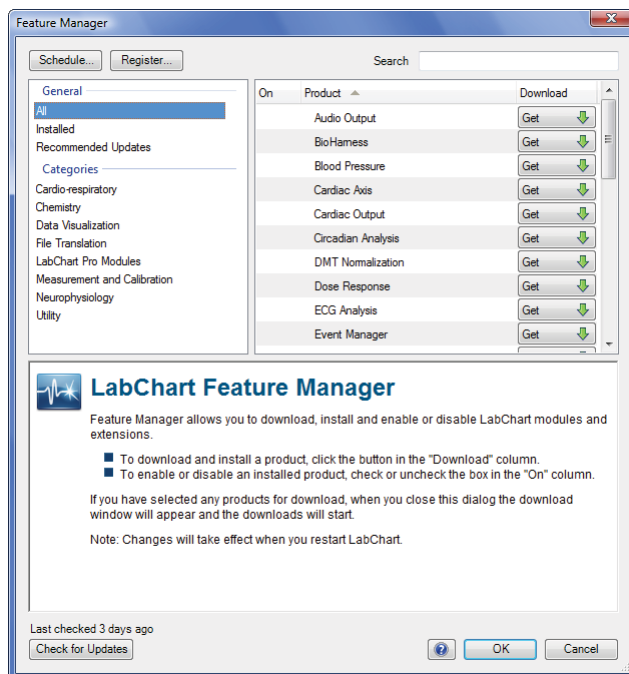

*Category (top left), Product (top right) and Information (bottom) Panels in Feature Manager Dialog*

## Navigating Around LabChart 7 Feature Manager

- Feature Manager is accessed by selecting **Help > Feature Manager**
- **Feature Manager** is divided into three panels:
  - **Product:** View the entire catalogue of LabChart software
  - **Category:** Group products from the Product panel into general and scientific application categories
  - **Information:** Describe the selected product major features and functions
- Find a suitable LabChart software by typing a keyword into the **Search** box
- Click the **Register** button to bring up the registration webpage
- **Check for Updates** to check for immediate software update checks or Schedule regular automatic software update checks


## Downloading and Installing LabChart Modules and Extensions

- In the **Product** panel, select a product by clicking on it. Product description will be displayed in the **Information** panel.

Product ▲	Download
Export MATLAB	Queued
Export Quick Time	Queued
Export WAV	Queued
Fast Response Output	Get
HRV	Get
Metabolic	Get
Multipoint Calibration	Get
Peak Analysis	Get
PVAN	Get

The 'Get' and 'Queue' button in the Product Panel

- Click the **Get** button to **Queue** for download. Repeat for other products.
- Click **OK**.
- Enter your registered user name and password in the **Connect to www.adinstruments.com** dialog.
- Click **OK**.

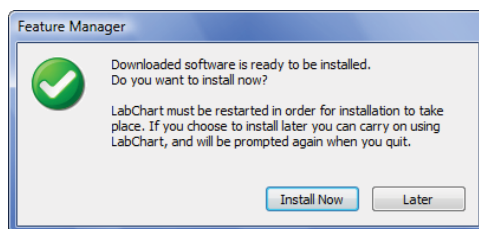

Option to either 'Install Now' or 'Later' downloaded software

- Choose to either **Install Now** software (LabChart would automatically shut down and restart) or **Later** to delay the installation process.


**Note:** A dialog would appear if an unlicensed LabChart Module is being installed. Click **Add License** to enter a purchased license code. Select **Buy Now** to purchase a license code or **Evaluate** to run the application in limited evaluation mode.

On	Product ▲	Download
<input checked="" type="checkbox"/>	Audio Output	
<input type="checkbox"/>	Dose Response	
<input type="checkbox"/>	ECG Analysis	
<input type="checkbox"/>	Spike Histogram	
<input type="checkbox"/>	Spirometry	
<input type="checkbox"/>	Video Capture	

Enabling and Disabling Installed LabChart software

- Select **Installed** from the **Feature Manager Category** panel. All installed products would be listed on the **Product** panel. Check/uncheck the box beside a product to enable/disable the product.

## Update LabChart Software

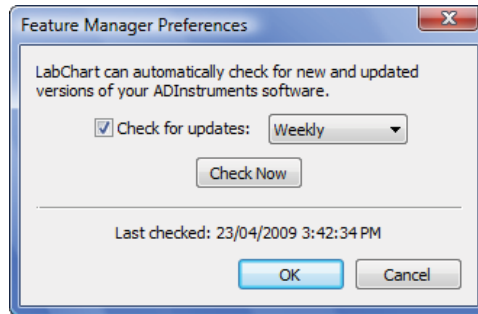

*Feature Manager Software Updates Dialog*

- Check for software updates from either **Help > Check for Updates**, **Check for Updates** button in **Feature Manager** Dialog or **Edit > Preferences > Feature Manager > Check Now**.
- Schedule automatic software update check from either **Edit > Preferences > Feature Manager** or **Schedule** button in **Feature Manager** Dialog.
- In **Feature Manager Category panel**, click on **Recommended Updates** to display updates available in the **Product** panel.
- Click the Update button on **Products** panel to **Queue** product for download.
- Click **OK**.